

SOCOREX®

ACURA® MANUAL PIPETTES




Fox
Scientific, Inc.

www.foxscientific.com • sales@foxscientific.com
(800) 369-5524



DWK LIFE SCIENCES
DURAN WHEATON KIMBLE

Acura® manual Pipettes

The Socorex® complete line of liquid handling products offers your lab a single source for superior quality pipettes, tips and accessories. Socorex pipettes which are manufactured in Switzerland are available exclusively by DWK Life Sciences (June 1st, 2017 WHEATON Inc was renamed to DWK Life Sciences) in the United States. Socorex *manual* pipettes meet and exceed users' needs for precision, accuracy and affordability. The Acura 825 and 835 are classic manual pipettes for all applications; the Acura 826 XS is the newest, light weight pipette designed intentionally to be shorter in length to facilitate working in small microtubes. In addition, all Acura *manual* pipettes are designed with optimal ergonomics to help prevent repetitive stress injuries. Specifically, the Acura 826 XS features an ultra-soft plunger which greatly reduces thumb and hand fatigue from long hours of pipetting.

Acura® manual 825

Product Advantages

- Smart and reliable volume adjustment
- Single-handed volume setting
- Precision digital display visible at all times
- Justip™ adjustable tip ejector* fitting most tips
- Swift-set user calibration system*
- Shock, UV-light and autoclaving resistance
- CE Certified IVD 98/79 EEC

* Socorex patented



825

- 0.1 - 2µL
- 0.5 - 10µL
- 1 - 10µL
- 2 - 20µL
- 5 - 50µL
- 10 - 100µL
- 20 - 200µL
- 100 - 1000µL




826 XS

- 0.1 - 2µL
- 0.5 - 10µL
- 1 - 10µL
- 2 - 20µL
- 5 - 50µL
- 10 - 100µL
- 20 - 200µL
- 100 - 1000µL

Acura® manual 826 XS

Product Advantages

- Excellent ergonomic shape and size ratio
- New springs and tightness seal for extra-smooth activation
- Reduced weight - lighter than leading brands
- Increased drivability with short, narrow shaft
- Easier to use when working in a fume hood
- Conical shaft end for easy reach in smaller microtubes

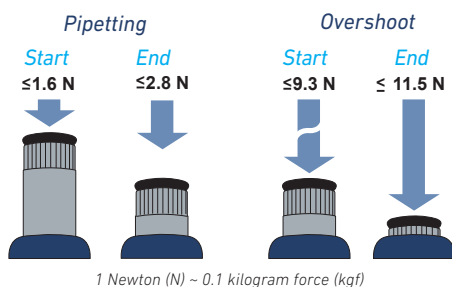

Extra Smooth Activation

- Unique tightness lip seal and new springs provide for ultra-soft pipetting, reducing hand fatigue while working.
- Forces measured on a 20 - 200µL model, are indicative of very limited finger efforts. Yet, the overshoot stop offers a clear tactile indicator.

Acura® manual Pipettes

Most Comfortable Pipetting 1

The Acura® manual fits everybody's hand with or without glove. Ergonomic shape, light weight and soft plunger travel care for better comfort.



Ejector Button 2

Ergonomically positioned, the large surface, soft padded ejector provides for low-pressure activation.



Adjustable Tip Ejection 3 4

Ejector shaft height adjustment system called Justip™ allows a wide selection of tips to tightly fit the nozzle.



Minimum Maintenance / Sterilization 5

Easy disassembling for thorough cleaning. Sterilization achieved by auto-claving instrument fully assembled (121°C / 250°F).



Instant Calibration 6 7

Swift-set user calibration system - with integrated key and locking mechanism - makes tedious procedures a thing of the past. Easy and precise, it is in full compliance with international standards and recommendations.



Safety Seal 8

Though easily removed when needed, the switch protection seal is autoclaving resistant.



NEW! Acura® manual 826 XS Pipette Starter Kit (W810200-XS)

Four pipettes, 4 racks of tips plus a stand

Special low price; Ideal for new labs!

(see back cover for a complete list of contents)

Calibration Seal Stickers

Cat. No.	Description	Qty/Case
W810620	Replacement seals	25
W810622	Replacement seals	100

Performance and Ordering Information - Acura® manual 825 and 826 XS Micropipettes

Cat. No.	Volume	Division	Inaccuracy (E%)			Imprecision (CV%)			Weight (g)	Tip Style
			Min. Vol.	Mid. Vol.	Max. Vol.	Min. Vol.	Mid. Vol.	Max. Vol.		
Acura 825										
W810300	0.1 - 2µL	0.002µL	< ± 6.0 % ¹⁾	< ± 4.0 %	< ± 2.0 %	< 5.0 % ¹⁾	< 3.3 %	< 1.5 %	88	10µL
W810302	0.5 - 10µL	0.01µL	< ± 2.5 % ²⁾	< ± 1.8 %	< ± 1.0 %	< 1.8 % ²⁾	< 1.2 %	< 0.5 %	88	10µL
W810304	1 - 10µL	0.01µL	< ± 2.5 %	< ± 1.8 %	< ± 1.0 %	< 2.5 %	< 1.6 %	< 0.7 %	88	200µL
W810306	2 - 20µL	0.02µL	< ± 2.5 %	< ± 1.8 %	< ± 1.0 %	< 1.7 %	< 1.1 %	< 0.5 %	88	200µL
W810308	5 - 50µL	0.1µL	< ± 1.5 %	< ± 1.3 %	< ± 1.0 %	< 1.0 %	< 0.7 %	< 0.4 %	90	200µL
W810310	10 - 100µL	0.1µL	< ± 1.5 %	< ± 1.2 %	< ± 0.8 %	< 1.0 %	< 0.6 %	< 0.2 %	92	200µL
W810312	20 - 200µL	0.2µL	< ± 1.5 %	< ± 1.1 %	< ± 0.6 %	< 0.6 %	< 0.4 %	< 0.2 %	90	200µL
W810314	100 - 1000µL	1µL	< ± 1.5 %	< ± 1.0 %	< ± 0.5 %	< 0.5 %	< 0.4 %	< 0.2 %	95	1000µL
Acura 826 XS										
W810300-XS	0.1 - 2µL	0.002µL	< ± 6.0 % ¹⁾	< ± 4.0 %	< ± 2.0 %	< 5.0 % ¹⁾	< 3.3 %	< 1.5 %	84	10µL
W810302-XS	0.5 - 10µL	0.01µL	< ± 2.5 % ²⁾	< ± 1.8 %	< ± 1.0 %	< 1.8 % ²⁾	< 1.2 %	< 0.5 %	84	10µL
W810304-XS	1 - 10µL	0.01µL	< ± 2.5 %	< ± 1.8 %	< ± 1.0 %	< 2.5 %	< 1.6 %	< 0.7 %	85	200µL
W810306-XS	2 - 20µL	0.02µL	< ± 2.5 %	< ± 1.8 %	< ± 1.0 %	< 1.7 %	< 1.1 %	< 0.5 %	85	200µL
W810308-XS	5 - 50µL	0.1µL	< ± 1.5 %	< ± 1.3 %	< ± 1.0 %	< 1.0 %	< 0.7 %	< 0.4 %	86	200µL
W810310-XS	10 - 100µL	0.1µL	< ± 1.5 %	< ± 1.2 %	< ± 0.8 %	< 1.0 %	< 0.6 %	< 0.2 %	87	200µL
W810312-XS	20 - 200µL	0.2µL	< ± 1.5 %	< ± 1.1 %	< ± 0.6 %	< 0.6 %	< 0.4 %	< 0.2 %	86	200µL
W810314-XS	100 - 1000µL	1µL	< ± 1.5 %	< ± 1.0 %	< ± 0.5 %	< 0.5 %	< 0.4 %	< 0.2 %	88	1000µL

Performance values obtained with double-distilled water at constant temperature (± 0.5°C) comprised between 20°C and 25°C in accordance with ISO 8655.

¹⁾ measured at 0.5µL ²⁾ measured at 1µL

Acura® manual Pipettes

Acura® manual 835 Macropipettes

Instruments combining high-tech materials with user-friendly ergonomics. Optimal for indoor and outdoor environmental analysis, clinical chemistry and cell culture applications, they offer superior performance and make pipetting safer than ever.


Product Advantages

- Smart and reliable volume adjustment
- Single-handed volume setting
- Interchangeable nozzle protection filter
- Justip™ adjustable tip ejector* fitting most tips
- Swift-set user calibration system*
- Shock, UV-light and autoclaving resistance
- Pasteur pipette fitting 2 and 5mL models

* Socorex patented



835
0.2 - 2mL
0.5 - 5mL
1 - 10mL

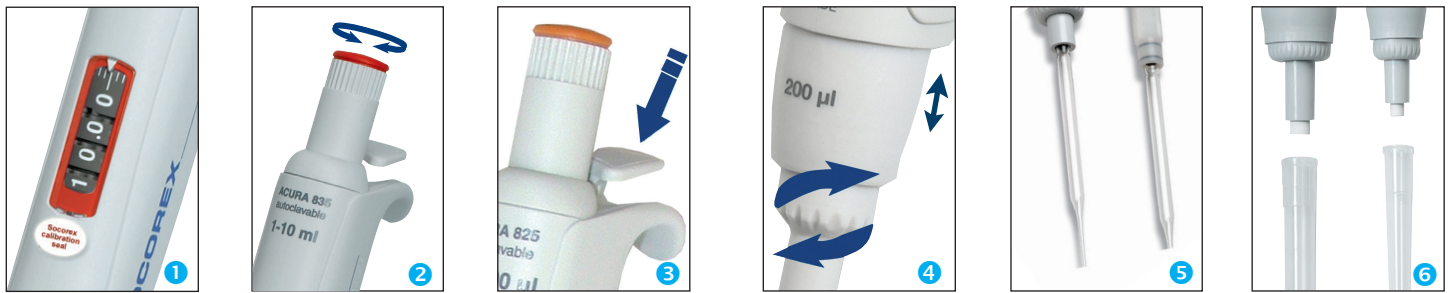

Display Always Visible 1

Located on the front of the instrument, the display window allows set volume to be visible at all times during pipetting. A comforting reassurance for the user.

Smart and Reliable Volume Adjustment 2

State-of-the-art micrometric volume setting is performed smoothly and precisely by turning push button. Finely machined click-stops and free rotating cap prevent any unwanted volume alteration.

Patented Tip Ejection System 3 4

Ergonomically positioned, the large surface, soft padded ejector provides for low-pressure activation. In addition, the shaft height adjustment system called Justip™ allows optimal distance setting between ejector and tip.

Pasteur Pipette Adapter 5

Adapter nozzles allow fitting of Pasteur pipette on Acura® manual 2 and 5mL models in addition to polypropylene tips. The adapters have two O-rings guaranteeing proper holding and tightness of standard 2mL Pasteur pipettes (Ø 7mm).

Nozzle Protection Filter 6

Protection against liquid overflow and contamination.

Performance and Ordering Information - Acura® manual 835 Macropipettes

Cat. No.	Volume	Division	Inaccuracy (E%)			Imprecision (CV%)			Tip Style
			Min. Vol.	Mid. Vol.	Max. Vol.	Min. Vol.	Mid. Vol.	Max. Vol.	
W810315	0.2 - 2mL	0.002mL	< ± 1.5 %	< ± 1.0 %	< ± 0.5 %	< 0.5 %	< 0.3 %	< 0.2 %	2mL
W810316	0.5 - 5mL	0.01mL	< ± 1.5 %	< ± 1.1 %	< ± 0.6 %	< 0.6 %	< 0.5 %	< 0.3 %	5mL
W810318	1 - 10mL	0.01mL	< ± 1.5 %	< ± 0.7 %	< ± 0.5 %	< 0.5 %	< 0.3 %	< 0.2 %	10mL

Performance values obtained with double-distilled water at constant temperature (± 0.5°C) comprised between 20°C and 25°C in accordance with ISO 8655.

Acura® manual Pipettes


Acura® manual 855 Multichannel Pipette

Built with the users in mind, these pipettes combine high-tech materials with exceptional, user-friendly ergonomics. They extend pipetting possibilities in 96-well microplates and guarantee outstanding performance and results.

Product Advantages

- 8 and 12-channel models up to 350µL
- Light weight and perfect hand fitting
- Precision digital display visible at all time
- Justip™ adjustable tip ejector* fitting most tips
- Swift-set user calibration system*
- 360° rotation of volumetric module

* Socorex patented

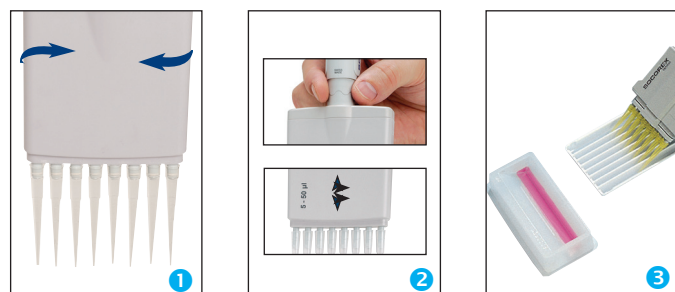

Optimal Working Position 1

Revolving over 360°, the volumetric module (lower assembly) helps reach the appropriate hand position.

Adjustable Tip Ejector 2

Position of tip ejector can be adjusted in no time within a 4mm span to perfectly match with tips used. Bowed shape of ejector head guarantees effortless, easy sequential tip removal.

Multichannel Reservoirs 3


Well adapted for multichannel pipettes, the reagent reservoirs offer various shapes and volumes.

8-Channel

- 0.5 - 10µL
- 5 - 50µL
- 20 - 200µL
- 40 - 350µL

12-Channel

- 0.5 - 10µL
- 5 - 50µL
- 20 - 200µL
- 40 - 350µL

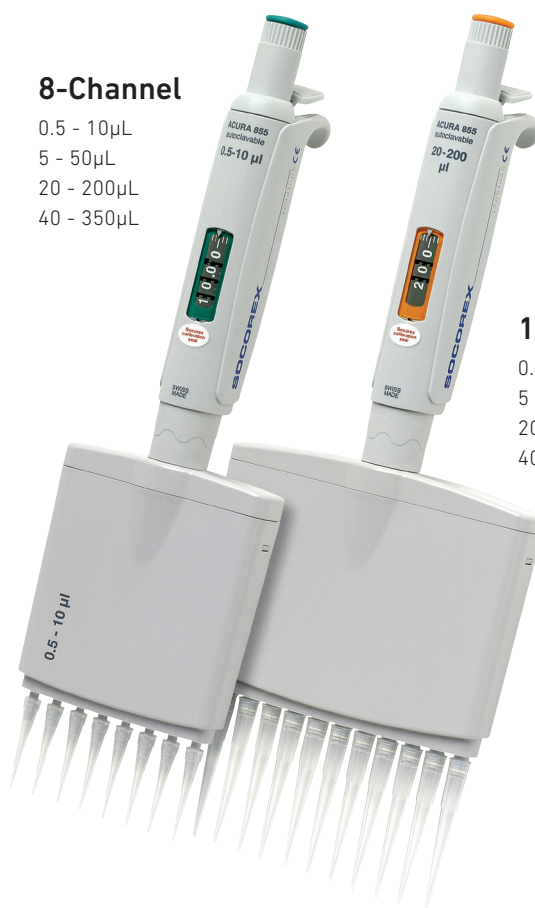

Performance and Ordering Information - Acura® manual 855 Multichannel Pipettes

Cat. No.	Volume	Division	Inaccuracy (E%)			Imprecision (CV%)			Weight (g)	Tip Style
			Min. Vol.	Mid. Vol.	Max. Vol.	Min. Vol.	Mid. Vol.	Max. Vol.		
8-Channel										
W810600	0.5 - 10µL	0.01µL	< ± 3.5 % ¹⁾	< ± 2.5 %	< ± 1.5 %	< 3.0 % ¹⁾	< 2.0 %	< 1.0 %	155	10µL
W810602	5 - 50µL	0.1µL	< ± 1.0 %	< ± 0.9 %	< ± 0.8 %	< 1.0 %	< 0.7 %	< 0.4 %	165	200µL
W810603	20 - 200µL	0.2µL	< ± 0.9 %	< ± 0.8 %	< ± 0.7 %	< 0.6 %	< 0.5 %	< 0.3 %	165	200µL
W810604	40 - 350µL	0.4µL	< ± 1.0 %	< ± 0.9 %	< ± 0.8 %	< 0.6 %	< 0.5 %	< 0.3 %	168	350µL
12-Channel										
W810606	0.5 - 10µL	0.01µL	< ± 3.5 % ¹⁾	< ± 2.5 %	< ± 1.5 %	< 3.0 % ¹⁾	< 2.0 %	< 1.0 %	184	10µL
W810608	5 - 50µL	0.1µL	< ± 1.0 %	< ± 0.9 %	< ± 0.8 %	< 1.0 %	< 0.7 %	< 0.4 %	199	200µL
W810609	20 - 200µL	0.2µL	< ± 0.9 %	< ± 0.8 %	< ± 0.7 %	< 0.6 %	< 0.5 %	< 0.3 %	201	200µL
W810610	40 - 350µL	0.4µL	< ± 1.0 %	< ± 0.9 %	< ± 0.8 %	< 0.6 %	< 0.5 %	< 0.3 %	203	350µL

Performance values obtained with double-distilled water at constant temperature (± 0.5°C) comprised between 20°C and 25°C in accordance with ISO 8655.

¹⁾ measured at 1µL

Acura® manual Pipette Accessories

Smartie Caps - Color Coding for Easy Identification



All Acura® pipettes have color-coded button caps for individual, lab or department identifications. Pick among 14 colors or choose the smartie mix for other color combinations.


Cat. No.	Color	Cat. No.	Color
Smarties Button Caps (6/pack)			
W820000	White	W820007	Rose
W820001	Grey	W820008	Pink
W820002	Vanilla	W820009	Purple
W820003	Lemon	W820010	Ice Blue
W820004	Yellow	W820011	Blue
W820005	Orange	W820012	Mint
W820006	Red	W820013	Green
Smartie Assortment (14/pack)			
W820014	Assorted colors (one of each above)		

Twister™ Pipette Stand, Model 336



Cat. No.	Color	Qty / Pack
W870152	Blue	1
W870154	Green	1
W870156	Orange	1
W870158	Grey	1
W870160	Red	1
W870162	White	1
W870164	Yellow	1

Nozzle Filters for Macropipettes

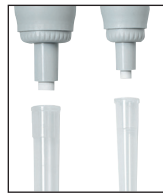

Add protection against liquid overflow or pipette contamination. Filters are interchangeable, non-sterile and non-autoclavable. Fits Socorex macro model pipettes.

Cat. No.	Fits	Material	Qty/Bag
W851342	Acura® manual 835, 2 and 5mL	PP fibers	250
851346	Acura® manual 835, 10mL	PP fibers	100

Work Stations, Models 337 & 340



337



340

Cat. No.	Color	Qty / Pack
Universal 7 Pipette Work Station, Model 337		
W851354	Ice blue	1
W851356	Mint green	1
W851359	Vanilla yellow	1
W851360	Pastel rose	1
W851361	Light grey	1
3 Pipette Work Station, Model 340		
W851369	Light grey	1

Pasteur Pipette Adapter



PVDF adapter fits pipette nozzle to accommodate standard glass Pasteur pipettes. The adapter has two FPM O-rings guaranteeing proper holding and tightness of standard glass 2mL Pasteur Pipettes.

Cat. No.	Description	Qty / Pack
W820125	For model 835, 2mL	1
W820023	For model 835, 5mL	1

Disposable Glass Pasteur Pipettes

Cat. No.	Length (mm)	Tip Length (mm)	Body Length (mm)	Shelf Pack	Qty/Case
357331	150	45	101	200	1000
357335	230	108	120	200	1000

Silicone Grease for manual Pipettes

Cat. No.	Description	Qty / Pack
W810624	1 gram Tube	5
W810626	10 gram Tube	1



Shelf Pipette Holder, Model 332





Cat. No.	Color	Qty / Pack
Shelf Pipette Holder, Model 332		
W844131	Transparent blue	2
W844132	Transparent blue	4



Qualitix® pipette tip selection



Ordering information


10 µL Tips	Size mm	Description		Sterile	Packaging		Cat. No.
		Colour	Style		Case	Quantity	
	31.4	Natural	10 µL ultra-microtips, fine pointed, graduated, autoclavable		1,000 / bag	1,000 tips	3090.0010B
		Natural	10 µL ultra-microtips, fine pointed, graduated, autoclavable		12x96 / rack	1,152 tips	3090.0010R
		Natural	10 µL <i>tipfill</i> ™ system incl. ultra-microtips, fine pointed, graduated		13x96 / case	1,248 tips	3090.0010T
				Empty rack for 10 / 200 µL tips to be used with <i>tipfill</i> ™ system, autoclavable	10 / case	10 racks	3070.0210ER
	71.0	Natural	10 µL capillary tips for gel loading, autoclavable		4x204 / rack	816 tips	302.0010GR


10 µL Filter Tips	Size mm	Description		Sterile	Packaging		Cat. No.
		Colour	Style		Case	Quantity	
	31.4	Natural	10 µL ultra-microtips, fine pointed, graduated	✓	12x96 / rack	1,152 tips	3090.0010FRS

10 µL Low Protein Binding Tips	Size mm	Description		Sterile	Packaging		Cat. No.
		Colour	Style		Case	Quantity	
	31.4	Natural	10 µL ultra-microtips, low binding, fine pointed, autoclavable		12x96 / rack	1,152 tips	3090.0010PR
		Natural	10 µL ultra-microtips, low binding, fine pointed, autoclavable	✓	12x96 / rack	1,152 tips	3090.0010PRS

20 - 100 µL Filter Tips	Size mm	Description		Sterile	Packaging		Cat. No.
		Colour	Style		Case	Quantity	
	49.6	Natural	20 µL filter tips, bevel point	✓	10x96 / rack	960 tips	3080.0020FRS
	49.6	Natural	100 µL filter tips, bevel point	✓	10x96 / rack	960 tips	3080.0100FRS

200 µL Tips	Size mm	Description		Sterile	Packaging		Cat. No.
		Colour	Style		Case	Quantity	
	49.6	Natural	200 µL universal tips, bevel point, graduated, autoclavable		1,000 / bag	1,000 tips	3070.0200B
		Natural	200 µL universal tips, bevel point, graduated, autoclavable		10x96 / rack	960 tips	3070.0200R
		Natural	200 µL <i>tipfill</i> ™ system incl. universal tips, bevel point, graduated		10x96 / case	960 tips	3070.0200T
	82.5	Natural	200 µL capillary tips for gel loading, autoclavable		6x96 / rack	576 tips	3080.0200GR
				Empty rack for 10 / 200 µL tips to be used with <i>tipfill</i> ™ system, autoclavable	10 / case	10 racks	3070.0210ER


200 µL Filter Tips	Size mm	Description		Sterile	Packaging		Cat. No.
		Colour	Style		Case	Quantity	
	51.2	Natural	200 µL universal tips, bevel point	✓	10x96 / rack	960 tips	3070.0200FRS


200 µL Low Protein Binding Tips	Size mm	Description		Sterile	Packaging		Cat. No.
		Colour	Style		Case	Quantity	
	49.6	Natural	200 µL universal tips, low binding, graduated, autoclavable		10x96 / rack	960 tips	3070.0200PR
		Natural	200 µL universal tips, low binding, graduated, autoclavable	✓	10x96 / rack	960 tips	3070.0200PRS


Qualitix® pipette tip selection


Ordering information



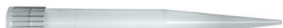
300 µL Filter Tips	Size mm	Description		Sterile	Packaging		Cat. No.
		Colour	Style		Case	Quantity	
	58.8	Natural	300 µL tips		10x96 / rack	960 tips	308.0300FR

350 µL Tips	Size mm	Description		Sterile	Packaging		Cat. No.
		Colour	Style		Case	Quantity	
	58.8	Natural	350 µL tips, autoclavable		1,000 / bag	1,000 tips	308.0350B
		Natural	350 µL tips, autoclavable		10x96 / rack	960 tips	308.0350R

1000 µL Tips	Size mm	Description		Sterile	Packaging		Cat. No.
		Colour	Style		Case	Quantity	
	88.9	Natural	1000 µL universal tips, fine pointed, autoclavable		1000 / bag	1,000 tips	3070.1000B
		Natural	1000 µL universal tips, fine pointed, autoclavable		6x96 / rack	576 tips	3070.1000R
		Natural	1000 µL <i>tipfill</i> ™ system incl. universal tips, fine pointed		5x96 / rack	480 tips	3070.1000T
				Empty rack for 1000 µL tips to be used with <i>tipfill</i> ™ system, autoclavable		6 / case	6 racks

1000 µL Filter Tips	Size mm	Description		Sterile	Packaging		Cat. No.
		Colour	Style		Case	Quantity	
	88.9	Natural	1000 µL universal tips, fine pointed	✓	6x96 / rack	576 tips	3070.1000FRS



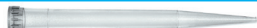






1000 µL Low Protein Binding Tips	Size mm	Description		Sterile	Packaging		Cat. No.
		Colour	Style		Case	Quantity	
	88.9	Natural	1000 µL universal tips, low binding, fine pointed, autoclavable		6x96 / rack	576 tips	3070.1000PR
		Natural	1000 µL ultra-microtips, fine pointed, autoclavable	✓	6x96 / rack	576 tips	3070.1000PRS

2, 5, 10 mL Macrotips	Size mm	Description		Sterile	Packaging		Cat. No.
		Colour	Style		Case	Quantity	
	115.5	Natural	2 mL macrotips, autoclavable		1x250 / bag	250 tips	851355
	123.2	Natural	5 mL macrotips, graduated, autoclavable		2x250 / bag	500 tips	W851362
		Natural	5 mL macrotips, graduated, autoclavable		2x50 / rack	100 tips	W851363
	150.0	Natural	10 mL macrotips, autoclavable		1x100 / bag	100 tips	851358
			Natural	Empty rack for 24x 10 ml macrotips, autoclavable		4 / case	4 racks

B = bag, E = empty, F = filter, G = gel, L = long, P = low binding, R = rack, S = sterile, T = *tipfill*™ refill system

Socorex® compatibility chart
















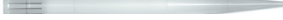

Chart indicates optimal combination of Qualitix® and Socorex® pipettes.


















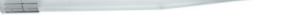

		Acura® manual and Acura® electro												Calibra® digital																		
		815					825 / 826 / 926					835 / 936		855 / 956		822	832	852														
		Up to 10 µL	10 µL / 15 / 20 µL	25 to 100 µL	120 to 200 µL	250 to 1000 µL	0.1 - 2 µL	0.5 - 10 µL	1 - 10 µL (V)	2 - 20 µL	5 - 50 µL	10 - 100 µL	20 - 200 µL	100 - 1000 µL	0.2 - 2 mL	0.5 - 5 mL	1 - 10 mL	0.5 - 10 µL	5 - 50 µL	10 - 100 µL	20 - 200 µL	40 - 350 µL	2 - 20 µL	10 - 100 µL	20 - 200 µL	100 - 1000 µL	0.2 - 2 mL	1 - 10 mL	1 - 10 µL	10 - 100 µL	20 - 200 µL	
10 µL Microtips	Tip ref.																															
		3090	•				•	•										•														
		302	• ²					• ²																								
	3090	•				•	•										•															
20 - 100 µL Microtips	Tip ref.																															
		3080	•					•	•	• ¹																						
	3080		•							•	•	• ¹																				
200 µL Microtips	Tip ref.																															
		3070	•	•	•			•	•	•	•	•							•	•	•											
		3080			•					•	•	•																				
	3070		•	•								•									•											
300 - 350 µL Microtips	Tip ref.																															
		308			•						•	•											•	•						•		
	308			•								•											•							•		
1000 µL Microtips	Tip ref.																															
		3070				•								•																		
	3070				•									•																		
2, 5, 10 mL Macrotips	Tip ref.																															
		851355												•														•				
		W851362													•																	
		W851363																														
	851358																															
	W835150																															

¹ Up to max. tip volume

Other brands compatibility

The Qualitix® line also fits major micropipette brands available in the market.

		Brand Such as models: Transferpette S, Electro					Eppendorf Such as models: Research Plus / Pro					Gilson Such as models: Pipetman, Classic, G, L				
		2.5, 10 µL	20, 50, 100, 200 µL	250, 300 µL	1000 (1200) µL	2, 5, 10 mL	2, 10 µL	20, 50, 100, 200 µL	300 µL	1000 µL	5, 10 mL	2, 10 µL	20, 50, 100, 200 µL	250, 300 µL	1000 µL	2, 5, 10 mL
10 µL Microtips	Tip ref.															
	3090	•					•					•				
	302	•					•					•				
	3090	•					•					•				
20 - 100 µL Microtips	Tip ref.															
	3080							• ¹					• ¹			
	3080		• ¹					• ¹					• ¹			
200 µL Microtips	Tip ref.															
	3070		•					•					•			
	3080		•					•					•			
	3070		•					•					•			
300 - 350 µL Microtips	Tip ref.															
	308			•									•			
	308			•												
1000 µL Microtips	Tip ref.															
	3070				•				•					•		
	3070				•				•					•		
2, 5, 10 mL Macrotips	Tip ref.															
	851355										•					
	W851362															•
	W851363															•
	851358															
	W835150															

		Rainin Such as models: Pipettite SL, EDP3 Plus					Sartorius Such as models: Proline Plus, Tacta, Picus					Thermo Such as models: Digital, F1, F2, F3, Novus					HTL, Labnet, Corning Such as models: Discovery, BioPette					
		2, 10 µL	20, 100, 200 µL	300 µL	1000 µL	5, 10 mL	2, 10 µL	20, 50, 100, 200 µL	250, 300, 350 µL	1000 (1200) µL	2, 5, 10 mL	2, 10 µL	20, 30, 50, 100, 200 µL	300 µL	1000 µL	2, 5, 10 mL	2, 10 µL	20, 50, 100, 200 µL	300 µL	1000 µL	5, 10 mL	
10 µL Microtips	Tip ref.																					
	3090	•					•				•				•							
	302																					
	3020	•					•				•				•							
20 - 100 µL Microtips	Tip ref.																					
	3080		• ¹					• ¹				• ¹				• ¹						
	3080							• ¹				• ¹				• ¹						
20 - 100 µL Microtips	Tip ref.																					
	3080		• ¹					• ¹				• ¹				• ¹						
	3080							• ¹				• ¹				• ¹						
200 µL Microtips	Tip ref.																					
	3070		•					•				•				•						
	3080		•					•				•				•						
	3070		•					•				•				•						
300 - 350 µL Microtips	Tip ref.																					
	308				•			•	•			•										
	308											•										
1000 µL Microtips	Tip ref.																					
	3070				•				•				•						•			
	3070				•				•				•						•			
2, 5, 10 mL Macrotips	Tip ref.																					
	851355																					
	W851362																			•		
	W851363																			•		
	851358															•						
	W835150																					

¹ Up to maximal tip volume

Brand names belong to their respective owners.

Acura® manual Pipettes

Acura® manual 825 / 835 Tripack

- Three adjustable micropipettes
- Large choice of volume combinations up to 10mL
- Nine different packs



Including: Sampling of pipette tips, silicone grease, Q.C. certificates and operating instructions

Cat. No.	Contents (Micro / Macropipette Models)		
W810319	2µL (0.1 - 2)	10µL (0.5 - 10)	50µL (5 - 50)
W810320	2µL (0.1 - 2)	20µL (2 - 20)	200µL (20 - 200)
W810321	10YµL * (1 - 10)	50µL (5 - 50)	200µL (20 - 200)
W810322	10µL (0.5 - 10)	100µL (10 - 100)	1000µL (100 - 1000)
W810323	10YµL (1 - 10)	100µL (10 - 100)	1000µL (100 - 1000)
W810324	20µL (2 - 20)	200µL (20 - 200)	1000µL (100 - 1000)
W810325	50µL (5 - 50)	200µL (20 - 200)	1000µL (100 - 1000)
W810326	100µL (10 - 100)	1000µL (100 - 1000)	5mL (0.1 - 5)
W810327	1000µL (100 - 1000)	5mL (0.5 - 5)	10mL (1 - 10)

* This pipette must be used with a 200µL pipette tip

Acura® manual 826 XS Pipette Starter Kit



Cat. No.	Contents	Qty
W810200-XS	W810302-XS (0.5 - 10µL)	1
	W810310-XS (0.5 - 10µL)	1
	W810312-XS (0.5 - 10µL)	1
	W810314-XS (0.5 - 10µL)	1
	Twister Blue Pipette Stand	1
	Rack with 10µL tips	1
	Rack with 200µL tips	1
	Rack with 1000µL tips	1

Acura® manual 826 XS TwiXS Pack

- Two adjustable volume pipettes
- Free shelf pipette holder
- Volume combinations up to 1000µL
- Six different packs



Including: Shelf pipette holder, pipette tips samples, silicone grease, Q.C. certificates and operating instructions

Cat. No.	Contents (Micropipettes)	
W810305-TXS	2µL (0.1 - 2)	20µL (2 - 20)
W810307-TXS	10µL (0.5 - 10)	100µL (10 - 100)
W810313-TXS	20µL (2 - 20)	200µL (20 - 200)
W810309-TXS	50µL (5 - 50)	200µL (20 - 200)
W810311-TXS	100µL (10 - 100)	1000µL (100 - 1000)
W810315-TXS	200µL (20 - 200)	1000µL (100 - 1000)



Fox Scientific, Inc.

www.foxscientific.com • sales@foxscientific.com
(800) 369-5524



Excellence in your hands

Contact us at [DWK.com](https://www.dwk.com)

© 2022 DWK Life Sciences • The trademarks used are owned by DWK Life Sciences GmbH, DWK Life Sciences LLC or DWK Life Sciences Ltd.
Acura®, Socorex®, Qualitix®, Twister™, Justip™ are trademarks or registered trademarks of Socorex Isba S.A.